

AstraZeneca 



# Automated Reading of Agar Plates using AI

24-25<sup>th</sup> March 2026

Andrew Gravett

Principal Scientist, Microbiology, Global Product Development, Pharmaceutical Technology & Development, AstraZeneca.

Reliable environmental monitoring plate reading powered by A.I.


**APAS** PharmaQC

*Never review a negative plate again.*




# Introducing APAS

APAS (Automated Plate Assessment System) Independence, by Clever Culture Systems (Adelaide, Australia), is an automated plate reader that uses a camera system and AI/ML model to count and sort plates.

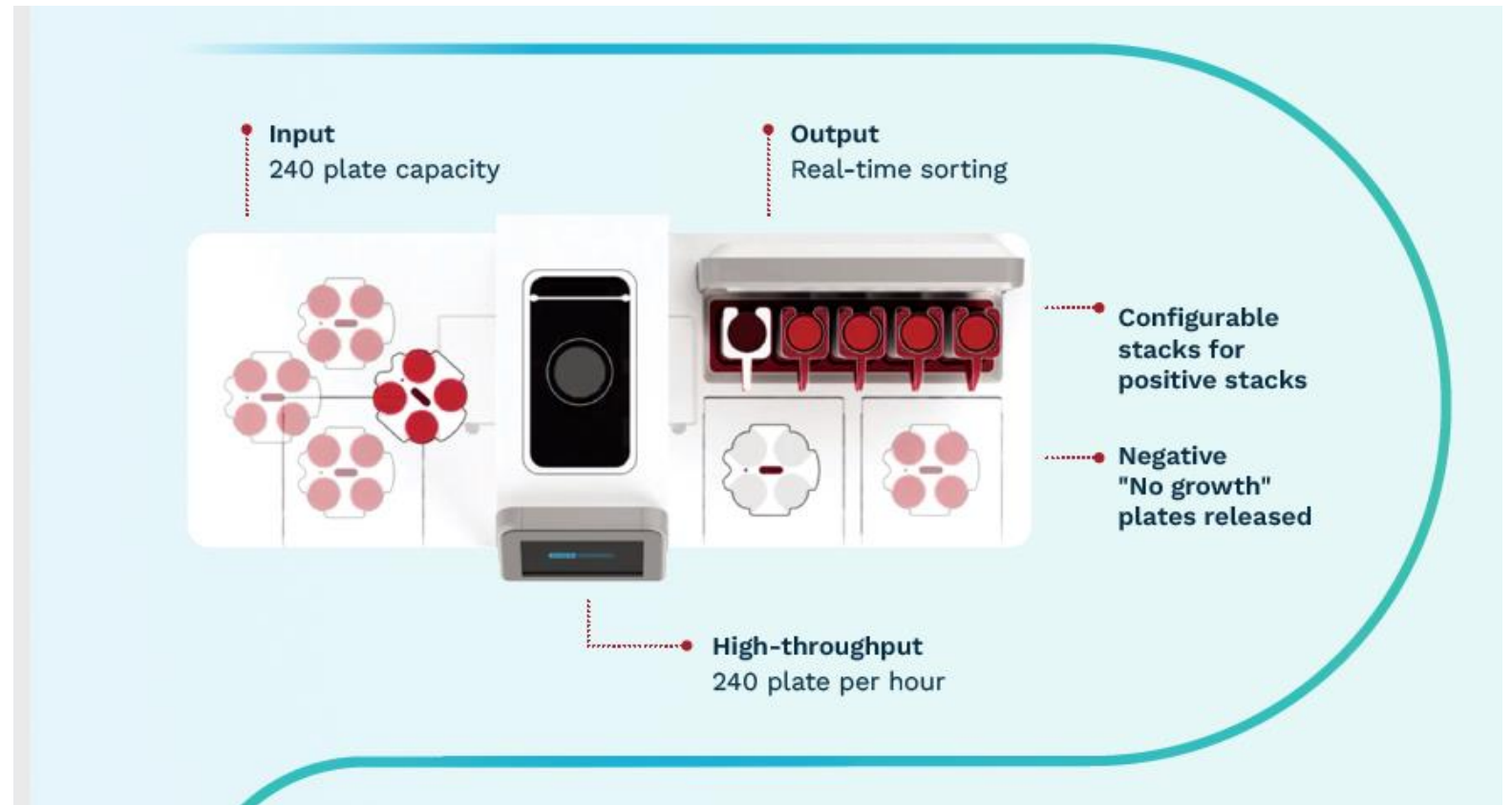

# Why were AstraZeneca interested?

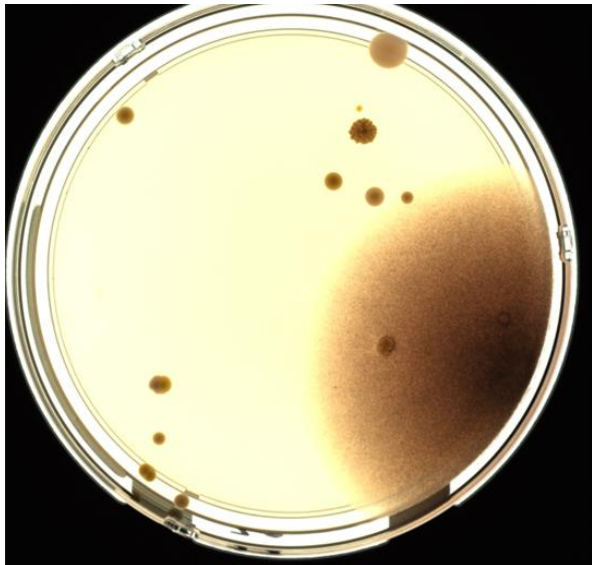

- Up to 30,000 agar plates are manually read and verified at large AZ sites
- Approximately 98% of these are negative (no growth)
- **A task requiring automation**
- Evaluated all available systems. Needed one which wouldn't require huge changes to existing infrastructure/processes:
  - Would work with numerous suppliers' petri dishes
  - Could cope with our numbers (without additional incubators)
  - Gave additional benefits of efficiency and data integrity
  - Could cope with different incubation protocols
  - Could save images/decisions if needed
- The solution was APAS – already approved use of AI/ML for clinical sample analysis.
- **We asked “Could AI/ML also be used to sort and count growth on Environmental Samples?”**



# How does APAS Work?




- **What is Artificial Intelligence in this instance?**
  - Basically a mathematical black box.
  - Algorithm developed to make decisions on images of agar plates
  - Based on the images it will decide if an EM agar plate is negative “No growth” or positive “Growth” or “contains artefacts”. Latter two sent for review by microbiologist.
- **But how does it decide?**
  - The black box is ‘trained’ (algorithm developed) to categorise what each pixel in the image is, based on a training set of images.
  - This is a long process of input pixel by pixel by microbiologists, and then testing the response to different ‘test’ plates/images.




# How is the algorithm developed?

- The algorithm is 'trained' in what the agar background looks like.
  - This is done by showing lots of images of different agar manufacturers agar plates and 'labelling' each artifact as background.
  - When APAS then detects these artefacts in real conditions, it will assign these pixels to the 'Agar Background' bucket.
  - The algorithm is then developed further by presenting it in training mode with lots of plates/images of different colony types.
  - It is then trained to put these colony types into a different segregation bucket nominally called 'White colony', 'Yellow colony', 'Filamentous Fungi', 'Transparent Colony' or 'Precipitate'
  - AZ shared images of 'real' environmental monitoring plates to train the algorithm in additional background and colony data.

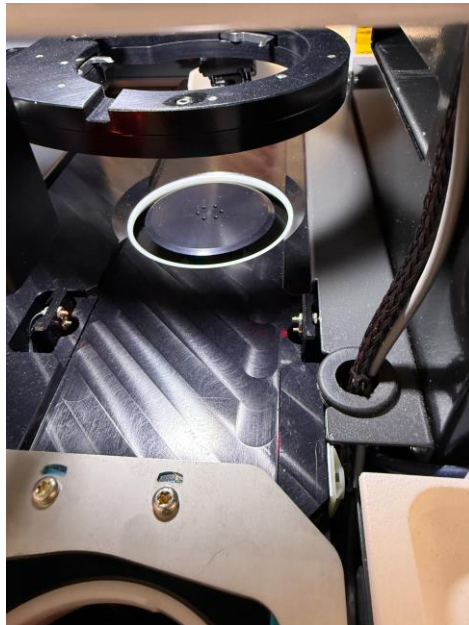

## Development of the Project and approach to validation and rollout

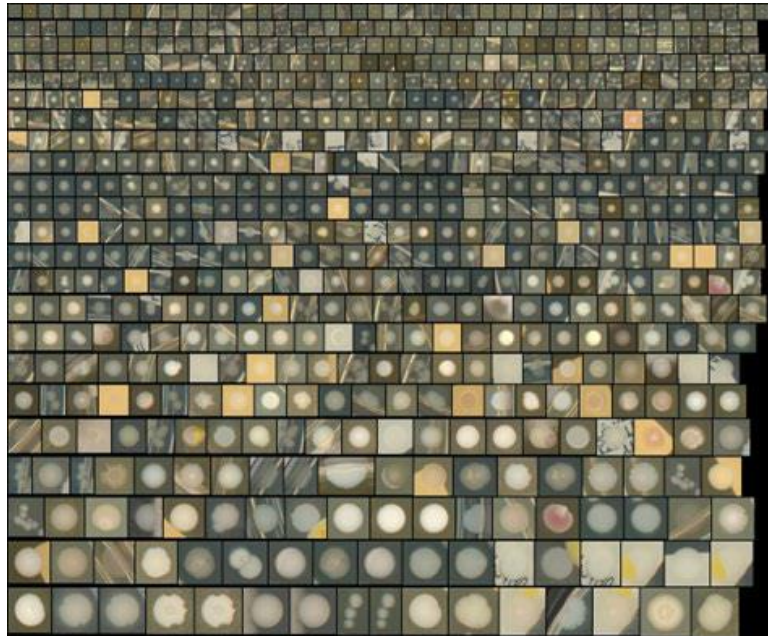

- Performed a proof of concept on an early prototype AI model
- Partnered with CCS to develop an Analysis Module to meet our automation needs of determining real microbial growth on agar plates and sorting them into separate stacks
- This can potentially save 80-90% of workload and bring other benefits such as:
  - Standardisation across global sites
  - Data integrity and data security
  - Ability to show images of actual plates to regulators
- Started with 90mm plates then repeated the process for 55mm plates once CCS developed the hardware
- CCS performed primary validation to USP<1223>, Ph Eur 5.1.6 and PDA Technical report No.33.
- AZ performed global OQ centrally and then rolled out to AZ sites around the world which then perform localised IQ and small PQ to save effort
- This process has taken 3 years but APAS is now in operation at 5 sites with 3 more at PQ and 2 more at IQ
- APAS is also now in use at other pharmaceutical companies and being evaluated by many more





# Obstacles

- Started the journey before AI was really known to the man on the street.....Would it be accepted?
- What is Raw Data in this case?
- Different views on Image Retention
- Different views on how to validate?
  - Equivalency versus target detection rate
- Lifecycle management
  - Need for periodic revalidation?
  - How to demonstrate no drift – ‘new isolates’
  - How to validate Analysis Module updates
- Have developed a Q and A guidance for common questions asked during interactions and conferences.
- Have Image retention policy.
- Have policy on what is considered raw data with APAS.



# What about Data Integrity?

- DI is a key benefit of automation
- Each agar plate has it's own individual barcode identifier
- Equipment software 21CFR part 11 compliant
- Validated automatic data transfer from APAS to LIMS system – removing error
- Images are retained for 45 days allowing review prior to batch release
- Much improved position compared to manual process which is documented and explained in a risk assessment




# Summary

- Long project with many hurdles
- Very worthwhile and more and more benefits and improvements are being seen following implementation
- Don't be afraid to innovate!

## 5 CLASSIFIER REVIEW - EXAMPLE IMAGES OF REMAINING ISSUES

The issues around the classifier can mostly be seen in these two examples:

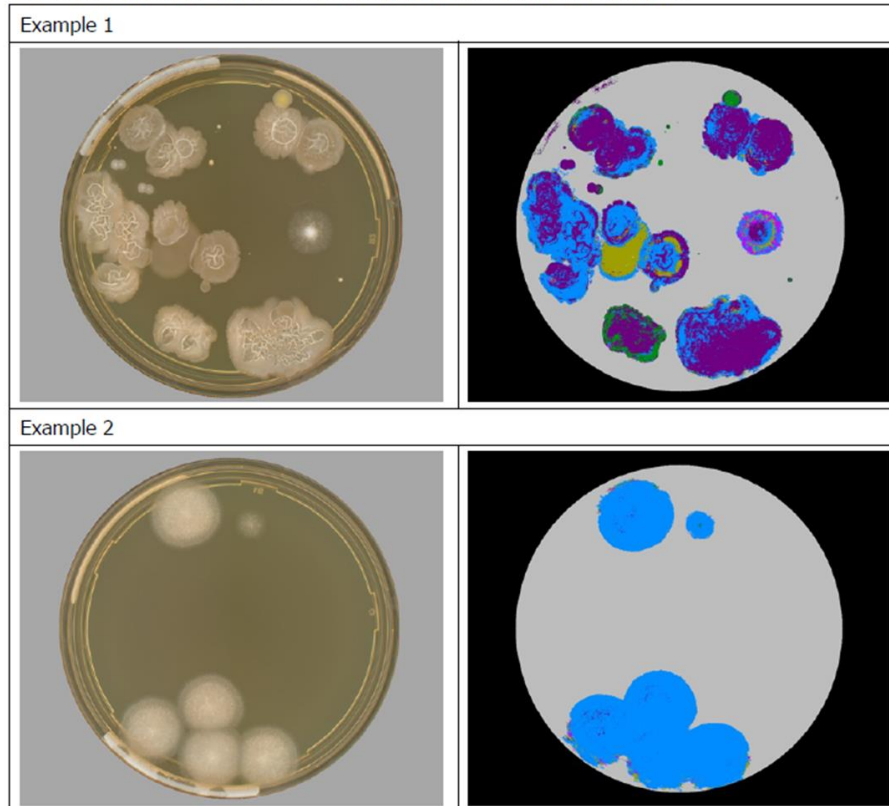

### Classifier 'Buckets'

- White colony
- Filamentous Fungi
- Yellow colony
- Transparent colony
- Precipitate

*Individual pixels are not enough to trigger a colony classification. A threshold value must be triggered for a particular 'bucket' to satisfy the algorithm that a colony is present.*



Thank you for your attention

**QUESTIONS ?**

### **Confidentiality Notice**

This file is private and may contain confidential and proprietary information. If you have received this file in error, please notify us and remove it from your system and note that you must not copy, distribute or take any action in reliance on it. Any unauthorized use or disclosure of the contents of this file is not permitted and may be unlawful. AstraZeneca PLC, 1 Francis Crick Avenue, Cambridge Biomedical Campus, Cambridge, CB2 0AA, UK, T: +44(0)203 749 5000, [www.astrazeneca.com](http://www.astrazeneca.com)

