

Mentype® AMLplus Profiler

Mentype® AMLplus Profiler is an analysis software program used in conjunction with the Mentype® AMLplus Mutation Analysis Kit from BIOTYPE GmbH. Data exported from GeneMapper™ 6.1 or GeneMapper™ ID-X 1.6 (Thermo Fisher Scientific) is converted by the Profiler into standardized PDF/TXT reports. The software runs exclusively locally on the laboratory's workstation; server operation or network access is not required.

Key Features at a Glance	
Single-user software with no server component	No application server, no database server, no service.
No network access required	The software does not communicate externally and does not require an internet connection. Neither a web server nor a network port is required.
No patient data	Processes only numerical fragment analysis data; no connection to personally identifiable information.
No need to exclude the software from virus scans	Common antivirus software generally does not cause conflicts.
Local admin privileges are sufficient	The installer writes to the program directory and the registry; no group-wide deployment or domain administrator is required.
No silent installation or GPO required	Standard installation via the Setup Wizard (GUI).
No online activation or licensing	The physical kits include a link and the password for downloading the installer. Internet access is required only to download the software; the installation itself runs in offline mode.
Clean uninstallation	No services or drivers are left behind when using the Windows "Programs and Features" function.

System Requirements (Verification Checklist) *Please check and tick off the following before installation*

- | | |
|---|--|
| ☐ Operating system: | Windows 11 (64-bit) |
| ☐ Prozessor: | Intel Core i5/ AMD Ryzen 5 or better |
| ☐ Memory (RAM): | At least 2 GB |
| ☐ Hard disk space: | At least 500 MB free (Installation + output files) |
| ☐ Screen resolution: | 1920 × 1080 or higher |
| ☐ PDF-Viewer: | Adobe Acrobat Reader or equivalent |
| ☐ GeneMapper™ 6.1/
GeneMapper™ ID-X 1.6 Software: | Already installed and running |
| ☐ Local admin rights: | Available for installation |
| ☐ Write permission: | User has write permissions for the desired output folder |

Note: Please verify system requirements against the current information in the product data sheet before installation

Data Access and Storage Locations	
Input:	GeneMapper™ 6.1 or GeneMapper™ ID-X 1.6 export file (TXT); manually selected by the user. The software does not have automatic folder access
Output:	Report files in PDF format and as a standardized TXT file for optimal reuse in LIMS systems. The user can freely choose the save location.
No writing outside the user folder	No system-wide write access during operation outside the installation directory.

If you have any technical questions regarding installation, our support team will be happy to assist you at support@biotype.de.