

Please follow the below guidelines to ensure the best performance



1. Purge from the thin needle at lower position

CAUTION: slow purging from **lower needle**, no more than **2 psi** purging for at least **10 minutes**. Argon (or other noble gas that is heavier than air) only.



2. Pull both needles out



3. Push the top cord gasket and apply clamp (make sure the **gasket is not bending when clamp is tight**)



4. Tight (keep squeezing the gasket via clamp, then tight it)

Apply clamp to push both ends of the cord gasket.

Follow these instructions will help maximize time of storing and safety.

Thank you for choosing PureAtom Scientific Container!

User Manual (in glove box, no purging noble gas needed)

Please follow the below guidelines to ensure the best performance



1. Pull both needles out after loading your sample, otherwise the container cannot be fully closed due to the pressure inside.



2. Push the top cord gasket and apply clamp (make sure the **gasket is not bending when clamp is tight**)



3. Tight (keep squeezing the gasket via clamp, then tight it)

Apply clamp to push both ends of the cord gasket.



Oxygen indicator will change the color to pink when O_2 level < 0.3% (left figure).

Follow these instructions will help maximize time of storing and safety.

Thank you for choosing PureAtom Scientific Container!