


## User Manual (PA-RT-002C)

Please follow the below guidelines to ensure the best performance:



1. Purge from the thin needle at lower position (skip this step when in glovebox)

CAUTION: slow purging from **lower needle**, no more than **2 psi** purging for at least **10 minutes**. Argon (or other noble gas that is heavier than air) only.

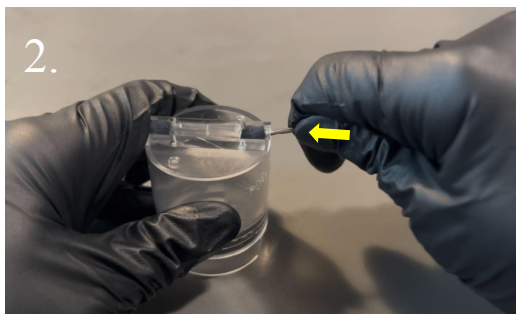

2. Pull both needles out and push in top gaskets to cover the holes



RT\_002C Video Guide

Scan the QR code to watch the video user guide:

We can also provide online demonstration. Please send your availability to [info.PA@pureatom.net](mailto:info.PA@pureatom.net) and we can answer your questions in the video call. Follow these instructions will help maximize time of storing and safety. Thank you for choosing PureAtom Scientific Container.

PureAtom LLC  
(208) 243-9610  
[info.PA@pureatom.net](mailto:info.PA@pureatom.net)  
6139 Bryson Lane  
Idaho Falls, ID 83402

