

User Manual

Please follow the below guidelines to ensure the best performance:




1. Purge from the thin needle at lower position (skip this step when in glovebox)

CAUTION: slow purging from **lower needle**, no more than **2 psi** purging for at least **10 minutes**. Argon (or other noble gas that is heavier than air) only.



2. Pull both needles out

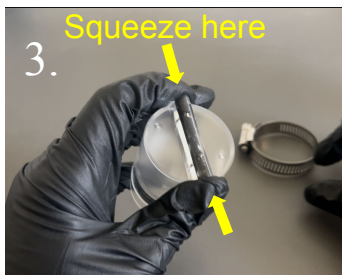

3. Push the top cord gasket and apply clamp (**make sure the gasket is not bending when clamp is tight**)



4. Tight (keep squeezing the gasket via clamp, then tight it)

Apply clamp to push both ends of the cord gasket.

Scan the QR code to watch the video user guide:



We can also provide online demonstration. Please send your availability to info.PA@pureatom.net and we can answer your questions in the video call. Follow these instructions will help maximize time of storing and safety. Thank you for choosing PureAtom Scientific Container.

PureAtom LLC
(208) 243-9610
info.PA@pureatom.net
6139 Bryson Lane
Idaho Falls, ID 83402

