

DissoPRO Software v3.6 Features

Pion's Rainbow R6 Concentration Monitor with in-situ fiber optic probes allows you to monitor dissolution rates in dissolution vessels without the need for sampling.

The fiber optic probes are submerged directly into the dissolution vessels and measure the UV absorption profile of your API as frequently as every 5 seconds and for as long as 28 days or more for extended release drug products.


The newly developed DissoPRO 3.6 software makes the integration into 21 CFR Part 11 - compliant environments straightforward.

Why Fiber optics?

In-situ real time fiber optics hold many advantages over traditional methods that involve sampling, fluidics/sample handling, and offline analysis.

In-situ Benefits:

- No filtration
- No pumping
- No sampling
- Dissolution profile with frequent data points (as quickly as every 5 seconds)
- Real-time dissolution curve for every channel, up to 8 channels

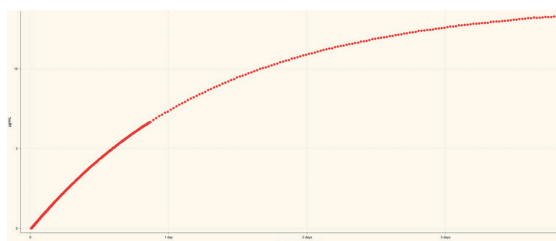


What does *in-situ* mean?

Fiber optic systems share a majority of elements with conventional UV spectrophotometers. The difference is that instead of drawing a sample and transferring into a cuvette and then the cuvette being inserted in the spectrophotometer for analysis, the *in-situ* fiber optic probes allow light to travel directly to and from the dissolution apparatus and analyses are performed directly in the dissolution baths within the windows of the fiber optic probes themselves. No sampling, no handling, no analysis delays.




Does it work in turbid media?


We provide a range of advanced mathematical approaches to correct the spectral data for turbidity including 4 different baseline correction methods and advanced second derivative spectrum analysis.

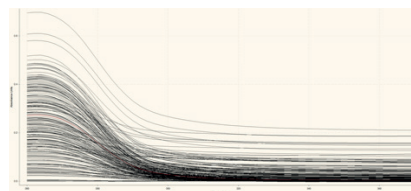

Direct spectrum; untreated

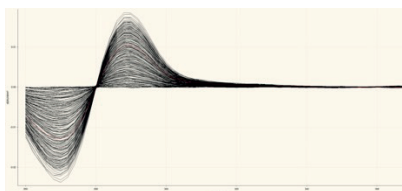


2nd derivative spectrum; corrected

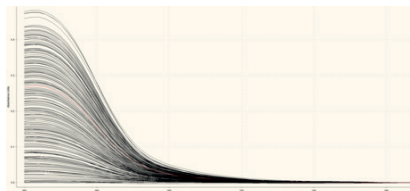


Direct spectrum; corrected

Software:

- Straightforward and powerful platform to collect dissolution data
- Can be integrated into 21 CFR Part 11 environment
- Key features of the software include: secure records, electronic signatures, audit trail, user security.

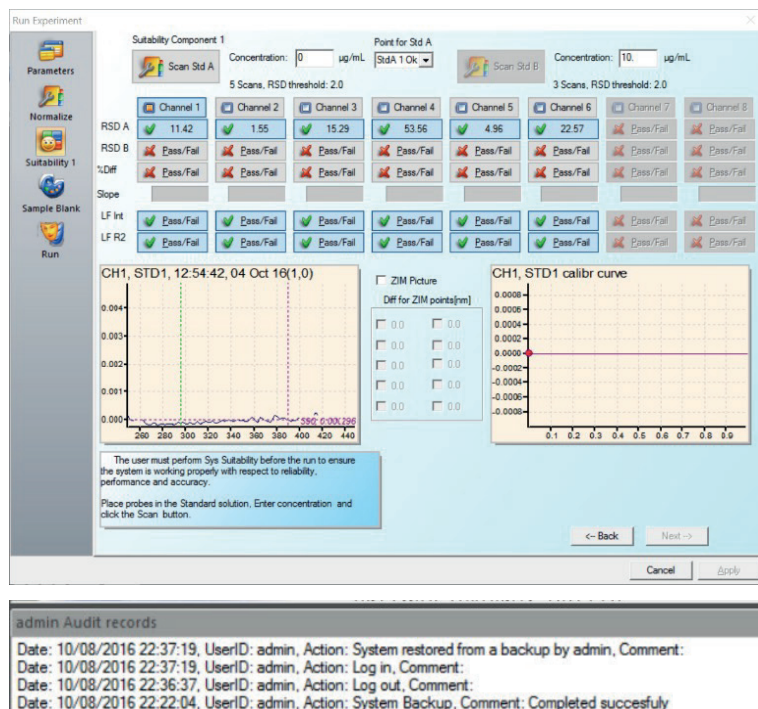


Field Title	Value
Project ID	TEST PROJECT
MethodID	000004
Method Title	Caffeine Demo 1
Author	dev1
CreationDate	17 May 2022 15:49:18
LastAccessDate	17 May 2022 15:49:18
Description	Caffeine test
Template Expirime	Empty_2022-05-17_15:39:32
Integrate Time Three	1000
hold	
Off threshold	50
Off threshold	FALSE
OPC Date	
Suitability Score A	5
Suitability Score B	5
Suitability Score C	5
Suitability RSD A	0.000000
Suitability RSD B	0.000000
Suitability RSD C	0.000000

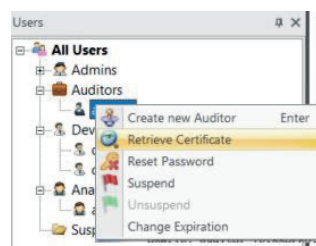


Integration - Working with your dissolution bath

- Compatible vessel covers are available (off-the-shelf or custom) whenever the bath design allows the integration
- Multiple probe designs are available (straight or J-style; fixed or variable pathlength) to optimize the data collection
- Broad pathlength range (1-20 mm) is available to fit the system to the desired concentration range

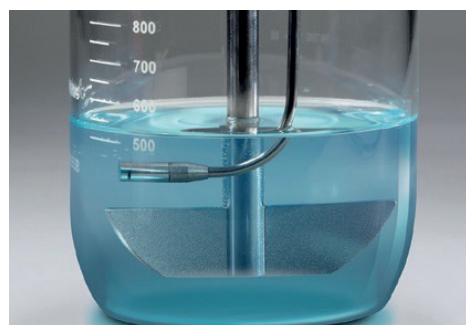

Documentation – fitting into your compliant environment

- Clear definition of features/functions relating to 21 CFR Part 11 are provided in the Design Specifications and Functional Specifications documents developed with a Computer System Validation specialist.
- Software and Hardware qualification is provided through detailed IQ/OQ/PV protocols.

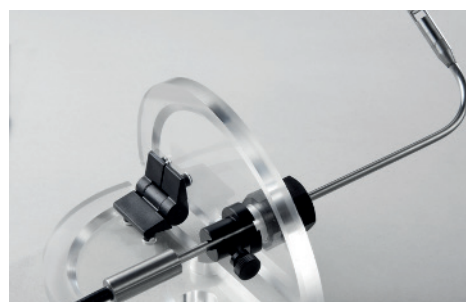

Pion stands behind the science

Pion Inc. | 10 Cook St. | Billerica, MA 01821 | MA | 01821 | +1 978.528.2020

Pion Inc. (UK) Ltd | Forest Row Business Park | Station Road, Forest Row | East Sussex RH18 5DW | +44 (0) 1342 820720

www.pion-inc.com | sales@pion-inc.com

An ISO9001:2015 company

Specifications subject to change without notice
Product sheet 90-40014