

SEP.T3.021 – Standard Procedure for the SiriusT3 Schedule to Load Manually


The procedure below will allow you to use the instrument without the Autoloader module:

- Please navigate to the following location on the instrument


PC: **C:\ProgramData\Pion\SiriusT3**

- If the Program Data folder is hidden:


- Navigate to “**Options**” in Windows Explorer

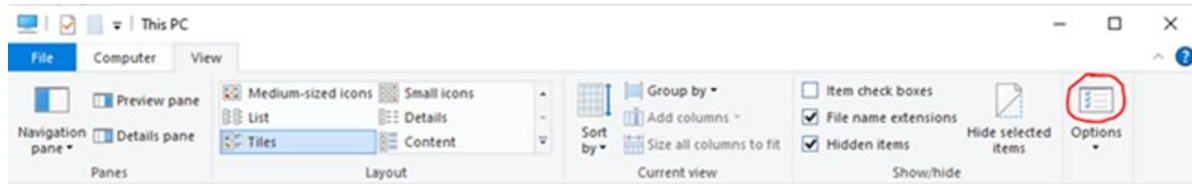


- In folder options they should scroll through the advanced settings of the “**View**” tab, locate “**Show hidden files, folders and drives**”, and then select the radio button for it.

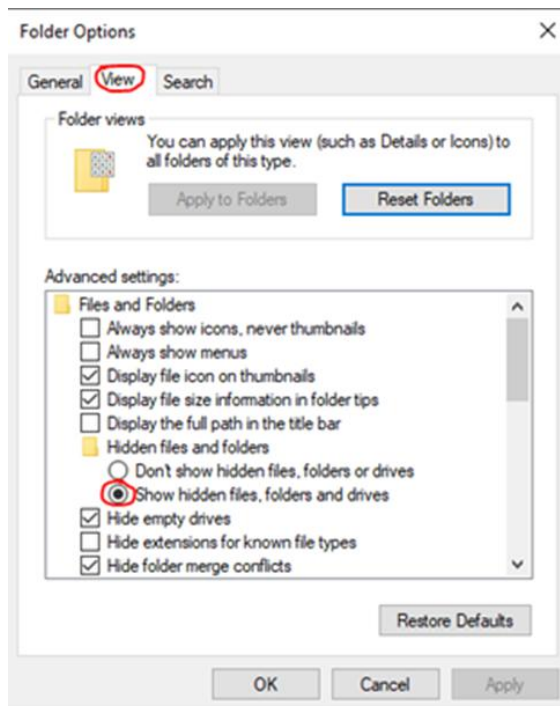

- Please take a copy of the **Instrument.XML** file and save it somewhere safe please.
- Open the **T3 Service** software and select the **Autoloader**.

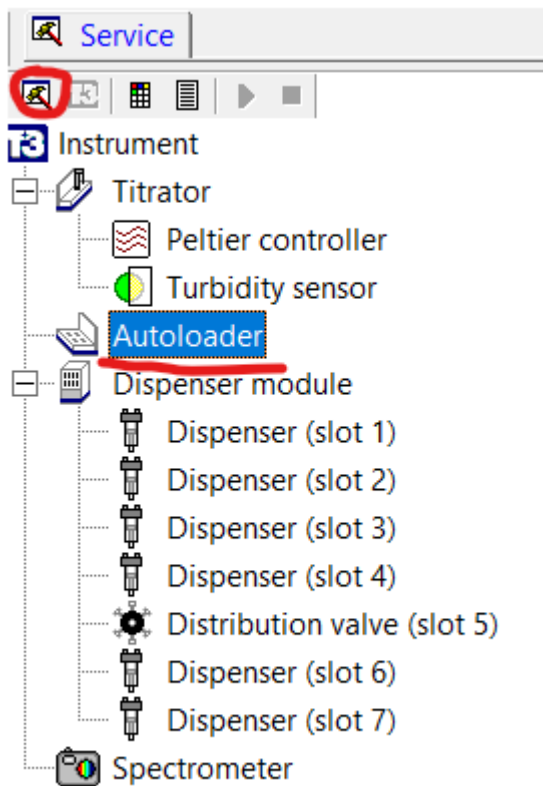

- With the **Autoloader** highlighted, press the **delete** button on the keyboard and click yes to Delete.
- Once deleted, please click the **Save** button. The Service text should now turn Blue.

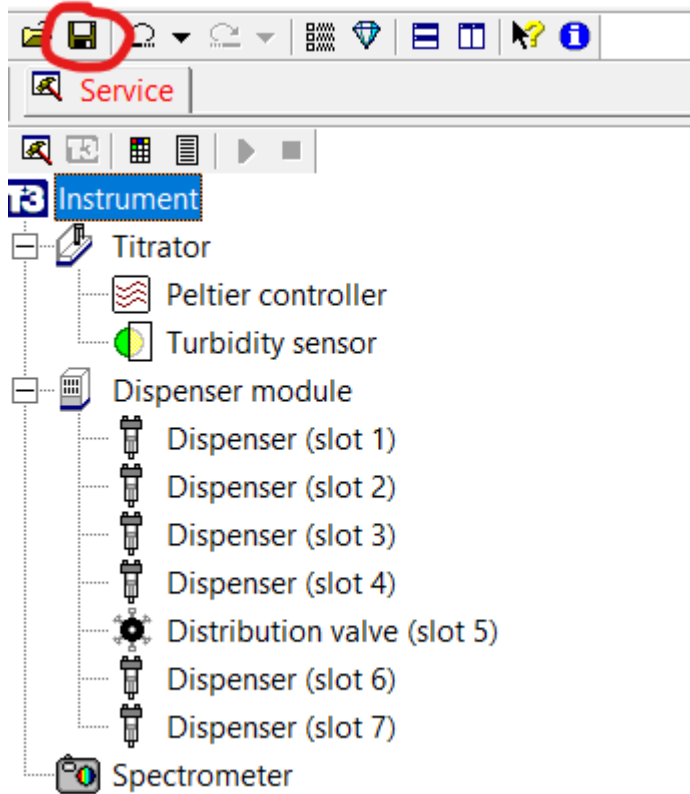

- Now when you open the Instrument software you can design the assays the same as before, but you won't have the Autoloader to run them. When you have designed the assay, you want to run, place an empty vial into the sample position and click the play button to run as normal. When the assay has finished remove the vial ready for the next assay.
- When the Autoloader is fully functional again, you can return the Instrument.XML file that you saved to a safe location (as this contains the Autoloader taught positions) to the Program Data location.