



Mach Zehnder Interferometer

Model: FOL - MZM


Interferometer is built by assembling the building blocks on a breadboard as per desired optical layout. Optics and laser head are held using kinematic mounts with two axis fine adjustment facility. Precision alignments for obtaining the interference fringes are done using kinematic mounts.

One of the simplest experiments using the interferometer is measurement of index of refraction of air with an airtight test cell placed in one of the optical arms of the interferometer. No other factor such as convective airflow, should be present, that can generate optical path differences.

The flexibility in fringe localization is one of the important advantages of the mach zehnder interferometer over other interferometers. Although many Mach-Zehnder interferometers use a rectangular arrangement, parallelogram arrangements are also possible. Vibration isolated supports for optical table is optional.

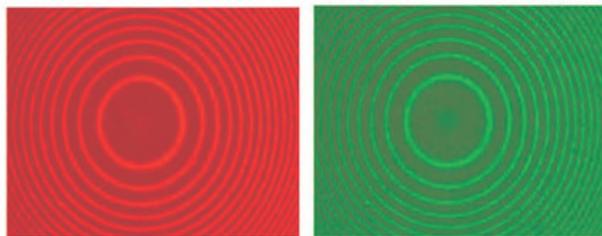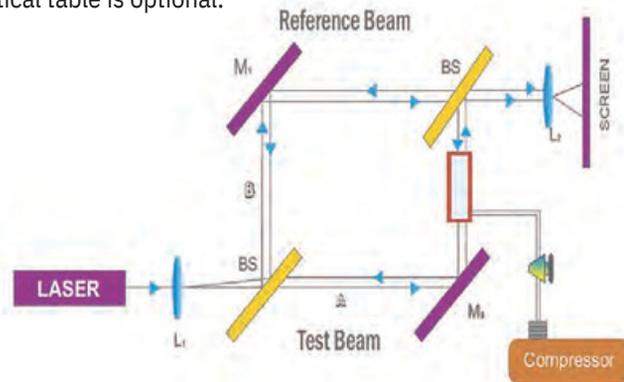

The Mach-Zehnder interferometer is a particularly simple device for demonstrating interference by division of amplitude. A light beam is first split into two parts by a beam splitter and then recombined by a second beam splitter. Depending on the relative phase acquired by the beam along the two paths the second beam splitter will reflect the beam with efficiency between 0 and 100%.

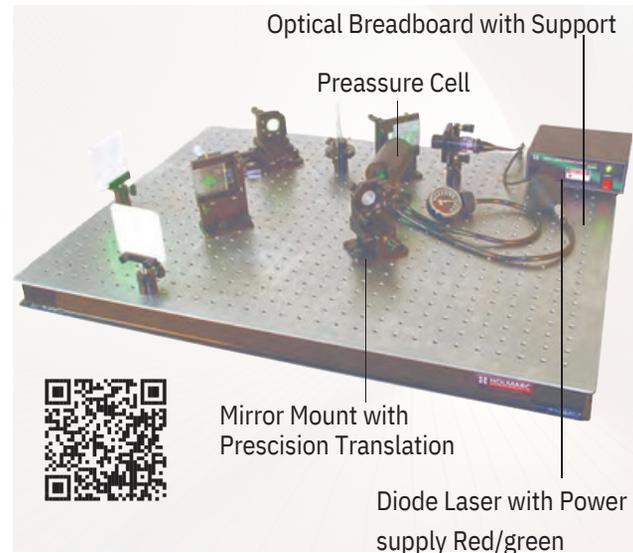

FRATURES

Kinematic mounts are used for fine movement of optical components

Diode laser is used as light source

All materials used in this setup are of laboratory grade with corrosion resistance

The optics used in this device are of research quality

Experiment Possible

To determine wavelength of laser beam.

The wavelength of laser is calculated by,

$$\lambda = (2d / N) \Delta$$

where 'd' is the change in position that occurs for 'N' fringes to pass and Δ is the calibration constant of the micrometer.

To find refractive index of a transparent material.

The light passes through a greater length of glass as the plate is rotated. The change in the path length of the light beam as the glass plate is rotated and relates the change in path length with the laser beam through air.

The refractive index of glass slide,

$$N = (2t - N\lambda) (1 - \cos\theta) / 2t (1 - \cos\theta) - N\lambda$$

Where t is the thickness of the glass slide, N is the number of fringes counted, λ is the wave length of light used and θ is the angle turned for N fringes.

To study refractive index change in air under different pressures and determine refractive index of air.

When a piece of material of thickness d is placed in one arm of the Michelson Interferometer, the change in optical path

length is given by $2dn$ where n is the difference in refractive index between the sample and the material it replaced (usually air). In other words, $2d(n - n_{air}) / \lambda$ extra wavelengths are introduced if air is replaced by a sample of refractive index n .

Scope of Supply

<p>Optical Breadboard with Support</p> <p>Model No. : ED-INT-12-BBS Dimensions : 800 mm x 600 mm Material : Stainless Steel Quantity : 1 No.</p>		<p>Kinematic Laser Mount</p> <p>Model No. : ED-INT-12-KLM Material : Black anodized Adjustments : Using 80 tpi lead screws Adjustment Range: +/- 3 degrees Quantity : 1 No.</p>	
<p>Beamsplitter Mount</p> <p>Model No. : ED-INT-12-BSM Degrees of freedom: 2 degrees Material : Black anodized Quantity : Aluminium alloy : 1 No</p>		<p>Mirror Mount with Translation</p> <p>Model No. : ED-IN T-12-MMP T Adjustment Range : +/- 3 degrees Material : Black anodized Least count of the micrometer : 0.01 mm Q uantity : 1 No.</p>	
<p>Mirror Mount with Precision Translation</p> <p>Model No. : ED-IN T-12-MMP T Adjustment Range : +/- 3 degrees Material : Black anodized Least count of the micrometer : 0.01 mm Q uantity : 1 No.</p>		<p>Rotation Stage</p> <p>Model No. : ED-IN T-12-R S Resolution : 2° / division Material : Black anodized Q uantity : Aluminium alloy : 1 No.</p>	
<p>Screen with Mount</p> <p>Model No. : ED-IN T-12-SM Dimension : 75mm x 75mm Q uantity : 1 No.</p>		<p>Pressure Cell</p> <p>Model No. : ED-IN T-12-P C Length : 10 cm Pressure Range : 0 - 300 mm Hg Q uantity : 1 No.</p>	
<p>Mirror with Cell</p> <p>Model No. : ED-IN T-12-MC Diameter : 25 mm Thickness : 6 mm Material : Borofloat Coating : Aluminium Q uantity : 2 Nos.</p>		<p>Beamsplitter</p> <p>Model No. : ED-IN T-12-B S Dimension : 50 mm x 50 mm Thickness : 6 mm R / T ratio : 50 / 50 Material : N-BK7 Coating : Aluminium Q uantity : 1 No.</p>	
<p>Glass Slide</p> <p>Model No. : ED-INT-12-GS Dimension : 75 mm x 25 mm Thickness : 1 mm Quantity : 5 Nos.</p>		<p>Diode Laser with Power supply (Green)</p> <p>Model No. : ED-INT-06-DLPG Wave length : 532 nm Optical power : 5 mW Quantity : 1 No.</p>	
<p>Diode Laser with Power supply (Red)</p> <p>Model No. : ED-IN T-12-D LPR Wave length : 650 nm Optical power : 5 mW Quantity : 1 No.</p>		<p>Diode Laser with Power supply (Red)</p> <p>Model No. : ED-IN T-12-D LPR Wave length : 650 nm Optical power : 5 mW Quantity : 1 No.</p>	