

INSTALLATION INSTRUCTIONS

REMOVAL OF THE PUMP MOTOR C-CLAMP

CAUTION: IMPORTANT STARTUP INSTRUCTIONS!

Tubing on all peristaltic pump assemblies on this machine have been protected during transportation and storage.

A metal C-clamp has been inserted to hold the spring rotor in each pump assembly. Please follow the steps below before powering up the machine.

1. Disassemble ALL pump assemblies from the machine by removing the holding screws.
2. Carefully remove the metal C-clamp from each clamp assembly.
3. Reconnect the Layall electrical connector.
4. Assemble the pump assembly back on the machine by replacing the removed holding screws.

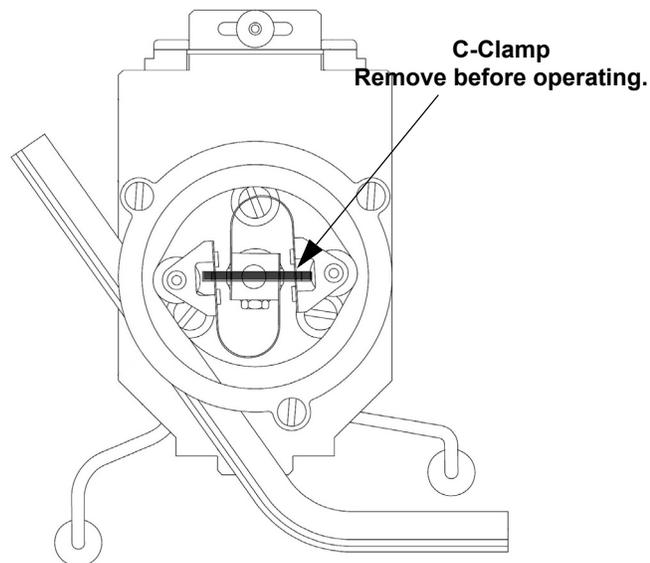


Figure 1