



When product integrity counts...

Manufacturers of laboratory calibrated temperature and humidity data loggers, real-time monitoring and cold chain solutions



Instructions for installation of TRW 6.3 software

- Admin rights are required to install software
 - Do not plug in reader until software is installed.
1. Go to Temprecord website: <https://temprecord.com/our-software/temprecord-windows-software/>
 2. Copy software from Downloads folder to PC
 3. Right click trw-set up exe and “extract all”
 4. Double click on set up file or EXE file and follow installation prompts.
 5. Plug in reader
 6. Re-booting may be required
 7. Open software
 8. Insert logger into reader (with logger facing up and hole exposed)
 9. Click space bar to access logger (The COM-PORT may need to be chosen)
 10. **Go to Parameters to enable use of the button on the logger**



For loggers that look like these, DO NOT use TRW software. For Mon-T2 data loggers please use [TempWrite software at:](https://temprecord.com/oursoftware/tempwrite/)

<https://temprecord.com/oursoftware/tempwrite/>